

Picking up of ultrathin section to grid is sometimes difficult. Collecting it by one hand, you may feel section escapes. Or you cannot get section where you would like to do. Especially when you use the grid with support film situation is much more complicated...

⇒ " MGH "solves your issues!!

How to use "Multi-Grid Holder?"




Put your grids into holders(up to 4 pieces).

Cut section is guided to 1st grid.

*Wherever you choose.

There is no specific rules.

Step 2
Collect 1st grid.

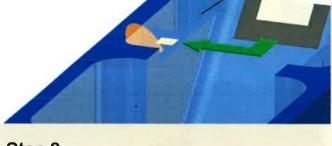

Step 3
Guide another section to 2nd grid.

Step 4

Collect 2nd grid.

*If you have 3rd (or also 4th) grid repeat above actions.

Description	Angle	width (mm)	Model No.	New	Exchange
Ultra (Wet)	35°	2.0 (2.0 to 2.2)	MGH 2035	OK	OK
		2.0 (2.0 to 2.2)	MGH 2045	OK	OK
*Recommended thickness range		2.5 (2.5 to 2.7)	MGH 2545	OK	OK
30 nm to 200 nm		3.0 (3.0 to 3.2)	MGH 3045	OK	OK

Remarks;

- 1. Replacement is only the same specification (especially diamond width) as yours.
- 2. For future improvement, some changes in appearance or specifications are done without notice.
- 3. For the preparation of very thin sections that are 30 to 40 nm, those results depend on your slicing circumstances.

*Such as sample width, embedding condition, hardness and so on.



All knives are made by...
SYNTEK CO., LTD.
Tel: +81(0)45-500-6603

Email: info@syntek.co.jp
Website: www.syntek.co.jp/en